

Description of options, upgrades and accessories for the laser beam stabilization system *“Compact”*

The basic configuration of the *Compact* laser beam stabilization system is fully equipped for stabilization tasks in laser machines or laboratory set-ups. It provides all required functions and parts (controller, detector/s, mirror holder/s and actuator/s, cables, etc.). However, for specific installations or lasers we offer a couple of adaptations or additional modules which further help to optimize the system for relevant parameters.

These adaptations and modules are described in this user manual. Some of these have to be integrated in the *Compact* system before the delivery. Others can be ordered as additional upgrades or accessories without the need for integration.

Contents

1. Controller modifications.....	2
1.1. Sample&hold circuit (“ADDA“) to fix the laser beam during laser off times.....	2
1.2. Adaptation for low repetition rates.....	2
2. Interfaces, additional inputs and outputs.....	3
2.1. External activation.....	3
2.2. Direct drive of Piezo actuators („Drive Actuator“.....	3
2.3. Voltage offset inputs to move the target position on PSDs (“Adjust-in“.....	4
2.4. Intensity outputs at controller (e.g. in case that UV or vacuum PSDs are used).....	4
2.5. Range outputs for monitoring applied Piezo voltages.....	5
2.6. Set-box for adjustment of target position on PSD.....	5
3. Detectors.....	5
3.1. Wide intensity detector: 4-quadrant diode with wide intensity range.....	5
3.2. 4-quadrant diodes for UV and IR.....	6
3.3. PSDs.....	6
4. Mirror mounts and mirror adapters.....	8
5. Vacuum adaptations.....	9
6. Optical components.....	9
6.1. Optical filters.....	9
6.2. Mirrors, lenses.....	10
7. Cables, other cable lengths.....	10
8. Shutter.....	10
9. Contact.....	11

1. Controller modifications

1.1. Sample&hold circuit (“ADDA“) to fix the laser beam during laser off times

In some applications with the laser beam stabilization the laser beam might be switched on and off during the operation. In laser off times there is no intensity on the detectors and hence no control signal for the closed-loop controller. In such situations, the *Compact* beam stabilization without additional measures will drive the Piezo-driven steering mirrors into a defined position, the so-called zero position (please refer to the *Compact* system's user manual for more information).

Once the laser is switched on again the stabilization will start its operation from this position. The zero position should have been used for the first adjustment of the optical set-up. That is why this is usually a good starting point for the stabilization. However, in cases of large drifts of the laser beam in the overall set-up, the zero position can – at least in the long term – strongly deviate from the required steering position. Switching on the laser after a time interval without laser beam and the resuming of the stabilization can therefore lead to an undesired initial spike of the beam position.

With the additional sample & hold circuit in the *Compact* beam stabilization system the positions of the steering mirrors can be fixed for an arbitrarily long time interval without control signal or laser intensity on the detectors. In that way it is possible to start the control-loop after switching on the laser not from the zero position but from that latest stabilized position.

The additional sample & hold (S&H) circuit is of special advantage for the following applications:

- In all systems where the laser must be switched on and off several times during the laser process, e.g. in material processing machines. Even if the system has drifted away from its basic adjustment, the beam position will start from the last stabilized position after resuming of the control loop. In this way an oscillation of the beam from the “zero position” to the desired position is eliminated and a potential faulty processing of the work piece is prohibited.
- In systems with a very large distance between the steering mirrors and the detectors. These set-ups bear the risk that a drift changes the adjustment in that way, that the laser beam will no longer hit the detector in uncontrolled intervals. Thus, it can happen, that the beam stabilization can not catch the beam on the detectors after a resuming of the stabilization when the laser was switched off for some time.
- In systems with very low repetition rates or lasers with irregular intervals of laser pulses (or pulse packages). If the S&H circuit of the beam stabilization is triggered for each laser pulse, the beam position will get closer to the desired position with each pulse.

The name “ADDA” is derived from the functional aspect that the actuators' drive signals are first AD converted and digitally stored before they are subsequently DA converted again and fed to the amplifiers of the mirror actuators. You can find more information about this module in a separate user manual.

1.2. Adaptation for low repetition rates

For lasers with low repetition rates (e.g. 10 ... 300 Hz) the detector electronics can be modified in order to maintain a position signal in the time gaps between two laser pulses.

Notes:

- Due to the low repetition rate the controller is then also optimized to low bandwidths.
- For lasers with even lower repetition rates we recommend our ADDA module (section 1.1).

2. Interfaces, additional inputs and outputs

Beside the inputs for the detectors and the outputs to the actuators the basic configuration of the *Compact* beam stabilization provides the following outputs:

- Status signal (see user manual of *Compact* system)
- Position signals x and y of each detector (analog voltage signal -5 to +5 V)

Other signal outputs or inputs can be provided as options.

Note: In some cases the arrangement of the connectors on the side panels have to be changed.

2.1. External activation

The external activation enables the change of the operation state of the beam stabilization system with an external signal.

There are three operation states. The specification of the control signal is as follows:

Signal (Level: 5V TTL)	Voltage range	Controller status	Reaction of <i>Active LED</i>
H (high)	2.4 – 5.0 V	Start	on
L (low)	0.0 – 0.8 V	Stop	off
Z (high impedance or not connected)		Manual mode according to selection on front panel	on/off

The external activation can be independently applied for stage 1 and stage 2 of the stabilization system. For this purpose, two LEMO connector plugs (series 00) are embedded on the right panel of the controller box. The inputs are marked as “Ext” and are next to the respective detector inputs.

2.2. Direct drive of Piezo actuators („Drive Actuator“)

As an option for the direct drive of the Piezo actuators (i.e. without feedback from the detectors) we can implement additional input channels to the controller. It is then possible to drive the actuators with an external voltage signal. This option makes use of the integrated 4-channel high-voltage amplifier of the system.

Specification:	Integrated 140V voltage converter
Inputs:	4 signal inputs (LEMO 00) on side panel, +/-5V
Outputs / to Piezo actuator:	4 voltage outputs on side panel, LEMO 0S series, 9V to 120V
Output impedance:	110 Ohm@1kHz, designed for high capacitive load
Power supply:	12V, 2A power supply

The input signal will be converted to a high-voltage signal which is fed to the Piezos.

Notes:

- The specification of the voltage range for the PKS / PSH Piezo actuators is -20V to +130V.

- The maximum tilt for the actuators is

Actuator type	Range
PKS	> 1 mrad (0.5 mrad in each direction)
PSH	> 2 mrad (1.0 mrad in each direction)
P4S30	> 4 mrad (2,0 mrad in each direction)

- We have specified the voltages for the valid range of the green range LEDs to values of 9V to 120V (max. range 0-130V).
- There is a non-linearity in both, the characteristics of the Piezos and the amplifiers. Therefore the signal will not be fully proportional to the input signal. If you need a precise and absolute position of the steering mirrors (without the control-loop which usually gives the position feedback) you should carry out a calibration of the angles versus voltages.
- It is also possible that the x and y axes of the same Piezo actuator vary strongly.

2.3. Voltage offset inputs to move the target position on PSDs (“Adjust-in”)

As described in section 3.3 the measurement principle of PSDs allows to move the target position on the detector by means of a voltage offset. For this purpose we can implement additional inputs for the x and y axes of both stages, 1 and 2. These inputs can be used to change the still stabilized beam position by an external source. The input voltage range of these inputs is -5V ... +5V.

Figure 1 shows the modified side panel of the controller box with additional “*Adj-in*” inputs for the voltage offset.



Figure 1: Left panel with four additional “*Adj-in*” inputs for x and y position of two PSDs

Note: The position vs. voltage characteristics of a PSD is usually not linear. Therefore, a calibration should be performed if the target shall be moved on a desired path.

2.4. Intensity outputs at controller (e.g. in case that UV or vacuum PSDs are used)

Since the UV and vacuum versions of the PSDs do not have the intensity level displays we can add additional intensity voltage outputs at the controller box. Figure 2 shows these outputs marked with “*Int*” close to the detector inputs. In this figure, the standard 4QD input connectors are labelled as input connectors for PSDs.



Figure 2: Right panel with PSD input connectors and intensity output connectors

2.5. Range outputs for monitoring applied Piezo voltages

In some applications it can be helpful to know the applied voltage ranges of the Piezos, e.g. to see whether or not the tilting range of the Piezos (and therefore the voltage range) is at its limits. If the Piezo actuators are combined with additional motorized mounts in order to enlarge the overall tilting range, the Piezo voltage can be used as a trigger to drive the motors.

2.6. Set-box for adjustment of target position on PSD

The set-box was developed as a simple tool for demonstration purposes and laboratory applications. It is an electronic box which offers potentiometers for x and y to manually adjust the target positions and outputs to the Piezo driver electronics (e.g. “Adjust In”, see section 2.3).

In an additional configuration it can be equipped with inputs for external signals so that an external signal from a signal generator can be added.

Note: The box can be equipped with x and y position and intensity displays. Please also ask for appropriate cables.

3. Detectors

3.1. Wide intensity detector: 4-quadrant diode with wide intensity range

In some applications the laser intensity is varied or modulated over wide ranges. The performance of the wide intensity detector is fully independent of the intensity. The signal amplification automatically adapts to changing powers. The power can vary by a factor of >1,000 without the need of exchanging the optical filters. External signals or user interactions are not required.

The detector has an excellent signal-to-noise ratio over the entire intensity range, so that the stabilization system always reaches the maximum resolution.

The function of the power level display is unchanged compared to the description in the *Compact* system's user manual, i.e. it can still be used to support the selection of the optical filters. In contrast, the potentiometer as described in section 4.5. of the user manual is omitted.

Note: Due to the wide intensity range it is possible to detect even lowest laser powers. Therefore, depending on the selection of the optical filters, the detection signal can be affected by ambient light.

3.2. 4-quadrant diodes for UV and IR

For lasers with UV and IR wavelengths we offer 4-quadrant-diodes with different sensitive areas and the following specs:

Specification:	UV 4-QD 3x3	IR 4-QD (InGaAs)	IR 4-QD (Ge)	Thermopile 4-QD
Wavelength range:	190-1,000 nm	900-1,700 nm	800-1,800 nm	190 nm-15 μ m
Sensitive area:	3x3 mm ²	$\varnothing = 3$ mm	$\varnothing = 5$ mm	18x18 mm ²

Figure 3 shows a photo of the sensor side of the UV 4-quadrant diode 3x3.



Figure 3: UV 4-quadrant diode with sensitive area of 3x3 mm²

Note: The Thermopile detectors are also suitable for CO₂ lasers at 10.6 μ m. Such lasers are often operated with powers in the range of several kW and large and heavy-weight mirrors. That is why the integration of the stabilization system can require additional measures for mounting the actuators and leaking a portion of the laser power to the detectors.

3.3. PSDs

As an alternative to our standard 4-quadrant diodes we offer PSDs (position sensitive device) for visible and for UV wavelengths with the following specs:

Specification:	VIS PSD	UV PSD
Wavelength range:	400-1,100 nm	200-1,100 nm
Sensitive area:	9x9 mm ²	10x10 mm ²

Compared to the 4-quadrant diodes the PSDs have a continuous measurement area. This leads to two possible advantages:

- 1) The sensitive area is not divided by a gap. Therefore, the PSD can be used in case of very small beam diameters or focused beams.
- 2) Whereas with the 4-quadrant diodes the target point is usually defined by their centre, in case of the PSDs any other point on the sensitive area can be chosen as a target point.

Applications:

If you use the PSDs instead of 4-quadrant diodes, the position detection is not limited to the centre as it is with 4-quadrant diodes. By adding a voltage to the signal of the PSD the target position where the laser beam shall hit the PSD can be moved. Still, the beam stabilization will provide full stabilization of the beam direction, but the direction itself can be manipulated. The external signal can be applied to the system via the “Adjust in” functionality described in section 2.3.

This feature can be used for different applications, e.g.:

- Place the PSDs before the laser is finally adjusted. Then adjust the laser and read-out the target position. Feed back the voltage for the new target position. The system will then stabilize the laser beam onto this position.
- Place the PSDs before the laser is finally adjusted. Then move the position on the detectors until you have the optimal laser adjustment.
- Move the laser beam to different points (or along a pattern) by moving the beam position on the PSDs. You can vary the laser beam direction with highest resolution and it is still stabilized.

Notes:

- A simple tool to move the target position by potentiometers is shown in section 2.6.
- The UV-PSDs do not have the position and intensity displays at the housing as the standard 4-quadrant diodes have.
- If we equip the beam stabilization with PSDs but no further measures, we use the electronic center (defined by a voltage of 0V for x and y position) as the target position.
- The position vs. voltage characteristics of a PSD is usually not linear. Therefore, a calibration should be performed if the target shall be moved on a desired path.
- The UV PSD can not be used for pulsed lasers with repetition rates below 1 kHz.

Please consider the additional input and output options of sections 2.3 and 2.4 which can further improve the functionality of stabilization systems with PSDs.

Adjustment of sensitivity

To make sure that the PSDs operate in the linear, controllable intensity range after a change of the laser intensity, you can adjust the signal gain by means of the digital potentiometers of each detector. In order to do this, please switch on the system (power ON) but do not activate the control loop (*Start/Stop* button switched off so that the green *Active*-LED and the LED besides the button do not shine). Then adjust the laser onto the detector and choose an appropriate gain with the digital potentiometer. The number of shining LEDs of the intensity display should be at least 3 and not more than 9.

The adjustment can be carried out with the delivered metal pin by means of soft pushing. There is a small push-button for each direction behind the bores in the housing (see the arrows in figure 4). With the upper push-button you can increase the gain step by step, with the lower push-button you can decrease it. There are 64 steps between the highest and the lowest gain. This corresponds to a change of the sensitivity by a factor of 20.

If you do not find a fitting adjustment, you can exchange the optical filters in front of the PSD sensors.

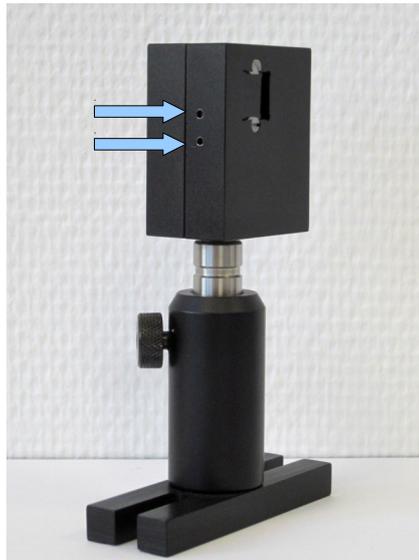


Figure 4: PSD detector. The arrows point onto the push-buttons of the digital potentiometers for the gain adjustment (which can be carried out with the delivered metal pin)

4. Mirror mounts and mirror adapters

Our standard mirror mounts have holders for 1" mirrors. The PKS mirror mount is also available for 0.5" mirrors.

Anyway, our beam stabilization system can also drive mirrors with larger dimensions. For such mirrors the system is equipped with specific Piezo-driven actuators and mounts, where the design is optimized with respect to control speed and tilting range. For 1.5", 2" and 3", the PSH and P4S30 steering mirror mounts can be equipped with appropriate adapters. Especially for mirrors with diameters of 2" or 3" we recommend the P4S30 mount. It works with four mutually clamped Piezo stacks and therefore yields a higher dynamic.

Figure 5 shows the P4S30 mount with two different mirror sizes. Figure 6 shows the PSH mount with a 2" mirror on the left side and a customised solution with a 4" mirror.



Figure 5: P4S30 mirror mount with a 1.5" mirror (left) and a 3" mirror (right)

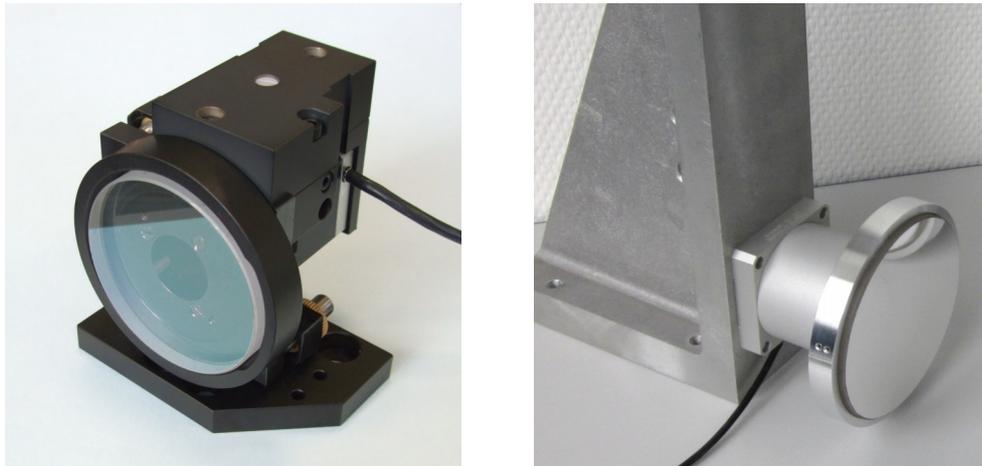


Figure 6: PSH mirror mount with a 2" mirror held by an adapter (left) and 4" mirror actuator (right)

Specification:	PKS	PSH	P4S30
Resonant frequency:	~ 700 Hz (1" mirror)	~ 840 Hz (1" mirror)	highest performance, even with large mirrors: ~ 1,200 Hz (1" mirror) ~ 400 Hz (1.5" mirror) ~ 300 Hz (2" mirror)
Mechanical tilt:	1 mrad (± 0.5 mrad)	2 mrad (± 1 mrad)	4 mrad (± 2 mrad)

Note: Please be aware that the bandwidths which can be achieved with larger mirrors is lower than with standard components.

5. Vacuum adaptations

Both, the detectors and the actuators can be adapted for use in vacuum. In case of the actuators, this is possible for vacuum pressures down to 10^{-11} mbar. But this is an extreme value. In case you intend to place some components in vacuum please let us know the conditions so that we can discuss and offer the required measures with you. Some measures (choice of materials and cables, sealing) are mainly focussed to avoid degassing and depend on the pressure. Other measures are important to protect the components themselves.

Note: The controller itself should not be placed in vacuum.

6. Optical components

6.1. Optical filters

We usually offer to integrate a pair of optical filters in front of each sensor. The filters have a size of 11.9×11.9 mm² and fit into the provided slot in the detector housing.

6.2. Mirrors, lenses

We can also offer appropriate filters or lenses for your set-up. However, since we do not produce optical components by our own we do not have any price advantages compared to our customers.

7. Cables, other cable lengths

The standard delivery of a *Compact* laser beam stabilization system includes all required cables to set up the system and to read out the positions. These are:

Cable set (included in standard delivery)	quantity	length
Detector → Controller	2	4 m (including the 3xMXC/power → Lemo adapter cable)
Actuator → Controller	2	PKS: 1.5 m (directly mounted to Piezo element)
		PSH: 1.2 m (directly mounted to Piezo element)
Actuator → Controller (Elongation)	1 pair for 1 actuator	10m
x-y-position cable (Lemo → BNC)	2	2 m

In addition to these cables we can also offer additional cables or cables with other lengths. The following table shows some examples.

Other available cables and/or lengths (examples)	Typical lengths
Elongation cables for 4QD detectors (Lemo → Lemo)	1 m ... 25 m
Elongation cables for actuators (Lemo → Lemo)	1 m ... 25 m
Cables for external activation (Lemo → BNC)	2 m
Cables to connect the intensity output (Lemo → BNC)	2 m

If you do not find the cable you need please do not hesitate to contact us. Since we assemble various cables in-house we can customize almost any cable and cable length.

8. Shutter

The laser shutter system *Beamblock* was especially developed for the use with our laser beam stabilization systems. It fulfils the safety function to block the laser beam in such cases where the stabilization system isn't capable to align the laser beam due to exceeding beam deviations. For this task the status signal "OK" of the stabilization system is used to control the laser shutter system externally.

The shutter system consists of a laser shutter and a shutter controller that allows to drive the shutter in three different modes of operation:

- confirm mode
- external mode
- manual mode

Thus the shutter can be controlled automatically by other devices or semi-manually with a safety confirmation by the user. Additionally it is possible for the user to open and close the shutter directly in the manual mode.

The input for external signals is prepared for standard TTL-levels, which allows the shutter system to be compatible to most other controllers (e.g. laser interlock or computer controlled devices).

Figure 7 shows photos of the laser shutter and the shutter controller.

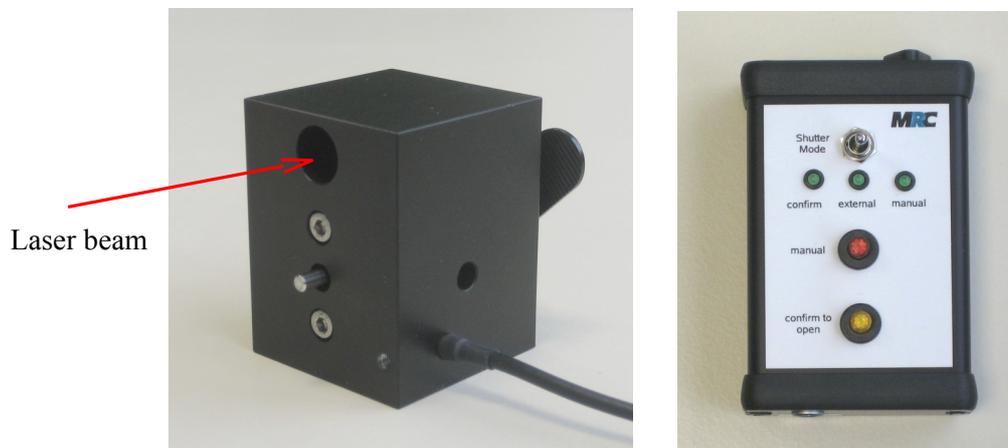


Figure 7: Laser shutter (left) and shutter controller (right)

You can find more information about this module in a separate user manual.

9. Contact

MRC Systems GmbH
Hans-Bunte-Str. 10
D-69123 Heidelberg
Germany

Phone: +49-(0)6221/13803-00
Fax: +49-(0)6221/13803-01
Website: www.mrc-systems.de
E-mail: info@mrc-systems.de